

Method

A variety of dairy products with a range of fat contents between 1% and 82% were analysed by NMR using an **MQC-23** fitted with a 26 mm hydrogen probe. Analyses were conducted using the **EasyCal** 'Fat in Foodstuffs' application software package.

Samples were prepared in duplicate by spreading 3–5 g of each sample on glass fibre filter paper and placing either a) on a metal dish or b) directly inside a glass vial. All samples were dried in an oven at 103°C for ~5 hours (drying time dependent on amount of sample). After drying, the open-filter samples were placed inside glass vials. All samples were conditioned at 40°C for one hour, prior to analysis. NMR measurement time was ~32 seconds.

The advantage of this method over extraction techniques is that a large number of samples can be prepared in batches.

Results

Figure 1 shows that a good calibration (Correlation Coefficient = 1.00; Standard Deviation = 0.57) can be obtained using all the samples of dairy produce and an empty tube which represents 0%.

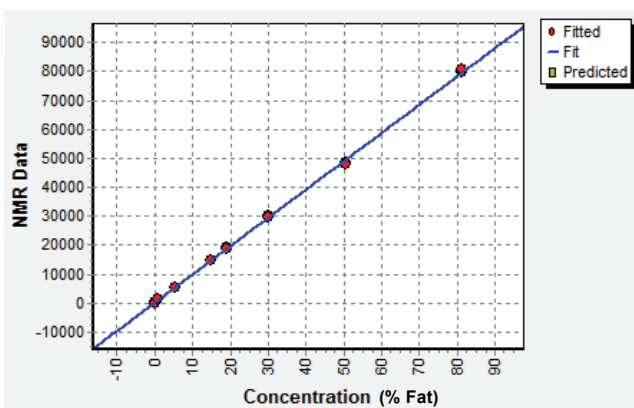


Figure 1

Since the hydrogen density of the fat components will vary with food type (e.g. butter, milk, vegetable oils and modified/ blended oils) it is recommended that a calibration is generated from fat derived from the same source material. Such a calibration may be generated by a 100% calibration method or a multi-point calibration covering the range of interest.

Sample	*Stated fat content (%)	Measured fat content (%)	Difference – relative to calibration line derived from all samples (%)
Butter	81.3	81.9	+0.6
Low fat strawberry yoghurt	0.8	1.1	+0.3
Raspberry yoghurt	5.5	5.3	-0.2
Low fat crème fraîche	15.0	14.9	-0.1
Crème fraîche	30.0	30.5	+0.5
Single cream	19.1	19.4	+0.3
Double cream	50.5	49.3	-1.3

Table 1: Fat content measured by NMR against the value *stated on the packaged produce.

Figure 2 shows an excellent correlation was achieved using four types of dairy produce and a zero control. By reducing the calibration range (e.g. 0-20%), data from the previous experiment gave a Correlation Coefficient of 1.00 and Standard Deviation of 0.23%.

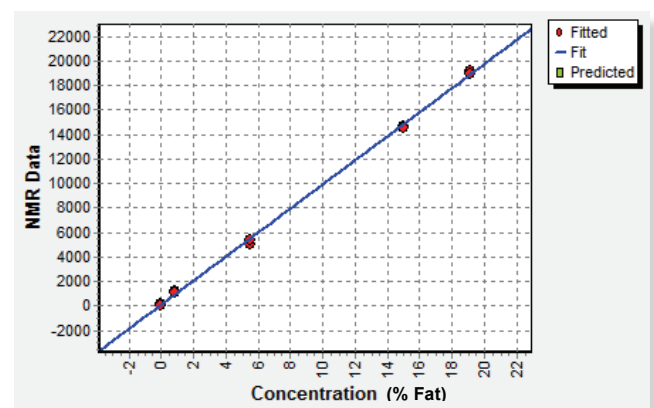


Figure 2



The Business of Science®



Conclusions

The **MQC-23** provides a means of rapidly measuring fat in dairy produce. Sample preparation is simple and conducive to very high sample throughput.



MQC – Enhanced Functionality

The instrument configuration used in this work may also be used to measure Oil in Dried Snack Foods, Fat in Chocolate and other Cocoa Derivatives. The functionality of the instrument may be increased to include the determination of Solid Fat Content (SFC), by the addition of a hydrogen probe (10 mm) and conditioning/tempering hardware as detailed in the AOCs/ISO SFC protocols.

Using Nuclear Magnetic Resonance to Test Fat Content in Foods

Case Study

How one contract laboratory dramatically increased sample throughput

The existing panel of business owners of Europe's leading independent testing laboratories is seeking alternatives to the standard solvent extraction methods. Established methods for testing the amount of fat in a range of foods. The contract laboratory with worldwide sites throughout the UK and Ireland, provides quality control and analysis services to the food industry. By converting from the wet chemistry method to **MQC** benchtop nuclear magnetic resonance (NMR) analysis for measuring fat content of foods, the lab reaped significant economic and environmental benefits. The box on the right is a brief overview of NMR's advantages over other standard methods.

Key Features of NMR

- Cost-effective: NMR's concentration range from 0.5 to 100 percent fat
- Precise calibration can be produced using a single fat sample
- Requires no sample preparation
- Sample measurement time is short
- Minimal sample preparation necessary
- No solvents are required
- Accurate for bulk measurements
- Accurate for sample granularity and product addition
- Non-destructive, facilitating repeatability measurements

Standard wet chemistry methods result in bottlenecks

Customer and product owners in the contract quality control sector are seeking alternative methods to the standard solvent extraction methods. Established methods for testing the amount of fat in a range of foods. The contract laboratory with worldwide sites throughout the UK and Ireland, provides quality control and analysis services to the food industry. By converting from the wet chemistry method to **MQC** benchtop nuclear magnetic resonance (NMR) analysis for measuring fat content of foods, the lab reaped significant economic and environmental benefits. The box on the right is a brief overview of NMR's advantages over other standard methods.

Oxford Instruments
The Business of Science®

Case Study: Fat in Food

Measurement of Oil Content in Dried Snack Foods

Application Note 2

Summary

- Up to 200 samples below ten min
- Chemically solvent free
- No hazardous waste produced
- Lowest most reliable technique available
- Suitable for difficult operators - simple, intuitive, visual software
- Simple linear calibration, no mathematics
- Suitable to sample from and addition
- Non-destructive, facilitating repeat

Advantages

- Up to 200 samples below ten min
- Chemically solvent free
- No hazardous waste produced
- Lowest most reliable technique available
- Suitable for difficult operators - simple, intuitive, visual software
- Simple linear calibration, no mathematics
- Suitable to sample from and addition
- Non-destructive, facilitating repeat

Advantages of NMR

- Cost-effective: NMR's concentration range from 0.5 to 100 percent fat
- Precise calibration can be produced using a single fat sample
- Requires no sample preparation
- Sample measurement time is short
- Minimal sample preparation necessary
- No solvents are required
- Accurate for bulk measurements
- Accurate for sample granularity and product addition
- Non-destructive, facilitating repeatability measurements

Oxford Instruments
The Business of Science®

No 2. Oil in Dried Snack Foods

Determination of Fat and Oil Content in Foodstuffs

Application Note 17

Advantages of benchtop NMR

- Cost-effective: NMR's concentration range from 0.5 to 100 percent fat
- Precise calibration can be produced using a single fat sample
- Requires no sample preparation
- Sample measurement time is short
- Minimal sample preparation necessary
- No solvents are required
- Accurate for bulk measurements
- Accurate for sample granularity and product addition
- Non-destructive, facilitating repeatability measurements

Oxford Instruments
The Business of Science®

No 17. Fat & Oil in Foodstuffs

Determination of Solid Fat Content in Edible Oils and Fats by the Official Direct Method (AOCS Cd 18b-93)

Application Note 13

Advantages of benchtop NMR

- Cost-effective: NMR's concentration range from 0.5 to 100 percent fat
- Precise calibration can be produced using a single fat sample
- Requires no sample preparation
- Sample measurement time is short
- Minimal sample preparation necessary
- No solvents are required
- Accurate for bulk measurements
- Accurate for sample granularity and product addition
- Non-destructive, facilitating repeatability measurements

Oxford Instruments
The Business of Science®

No 13. Solid Fat Content (SFC)

visit www.oxford-instruments.com for more information

www.oxford-instruments.com

This publication is the copyright of Oxford Instruments and provides outline information only which (unless agreed by the company in writing) may not be used, applied or reproduced for any purpose or form part of any order or contract or be regarded as a representation relating to the products or services concerned. Oxford Instruments' policy is one of continued improvement. The company reserves the right to alter, without notice, the specification, design or conditions of supply of any product or service. Oxford Instruments acknowledges all trademarks and registrations. Microsoft and Windows are registered trademarks of the Microsoft Corporation in the United States and other countries. Ref. LN_010_06_12



The Business of Science®

